


Guide for Identifying and Correcting Common Cell Growth Problems with Adherent Cells

CORNING

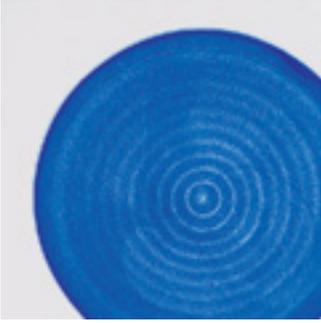

Guide for Identifying and Correcting Common Cell Growth Problems with Adherent Cells

Table of Contents

Introduction.....	3
Surface Treatment Process.....	3
Problems Related to Technique	4
Problems Related to Incubators	7
Problems Related to Culture Media	10
Problem Solving Suggestions.....	12
Acknowledgment	13
References	14

Acknowledgment

Corning wishes to acknowledge the contributions of John Ryan to this article.

Introduction

Although very popular among life science researchers, cell culture can be a very difficult tool to work with in the laboratory. Unlike other common laboratory tools such as electrophoresis or chromatography, cell culture makes dynamic, ongoing use of living organisms. These living cultures often respond to our mistakes not only by the erratic behavior common to other laboratory tools, but by dying – the total, irreplaceable destruction of the tool itself.

As one of the major suppliers of vessels, media, and sera used in cell culture, Corning Life Sciences often receives calls from anxious customers who are experiencing growth or attachment problems with their cell lines to determine what is responsible for the erratic behavior or loss of their cultures. Usually the customer, searching for a cause (and solution) to their problems, suspects that a change or mistake in the manufacturing process is responsible, or that there is an issue with the culture media or sera, or that perhaps the special surface treatment process used on most cell culture vessels was not properly done.

Because of the complex nature of cell culture, identifying the underlying causes of culture behavior problems is often a difficult, time-consuming task. Erratic culture behavior can take many forms; unusual growth patterns or inconsistent, spotty, and uneven cell attachment are the most common problems. Gradual or abrupt changes in growth rate or unexplainable experimental results are also experienced occasionally. In cell culture, any sudden change is suspect and a potential problem and, therefore, to be avoided.

Corning Life Sciences has spent many years helping customers deal with these problems. By using some of the information gained from these customer contacts, Corning has produced this guide to help identify and solve some of the common and uncommon causes of cell culture problems. The focus will be limited to three of the most common problem areas: technique, incubators, and media. In addition to the information provided in this guide, it is strongly suggested that you refer to the articles listed in the references or visit www.corning.com/lifesciences for additional help and recommendations.

Technical Assistance

For assistance in this area, please contact Corning Life Sciences Scientific Support at 800.492.1110, +1.978.442.2200, or email ScientificSupport@corning.com.

Surface Treatment Process

The initial and most common suspects for cell culture problems are usually either the culture vessels or the medium being used. Problems associated with culture medium will be dealt with in a later section. Much of the suspicion surrounding plastic cell culture vessels is due to the mystery or lack of understanding concerning the special treatment process used to modify the surface of the plastic. Virgin polystyrene, the resin used to manufacture most cell culture vessels, is hydrophobic in its untreated state. Proteins important for cell attachment, such as fibronectin and vitronectin (found in media containing serum) do not bind well to this natural surface resulting in poor cell adhesion and growth. Keep in mind, if you are going from serum containing culture media to serum-free media, this will impact the attachment of some cell types. Because of this, either a corona discharge or plasma treatment is used under carefully controlled conditions during the manufacturing process to insert oxygen atoms (in the form of carboxyl, hydroxyl, and ketone groups) shown in Figure 1, into the backbone of the polystyrene chain^{1,6,14}. This is the standard surface treatment used for most polystyrene culture vessels. This alteration of the plastic polymer (not a peelable coating) results in a hydrophilic surface with a net negative charge that creates a surface suitable for cell attachment and growth. The culture vessels are then sterilized and thoroughly evaluated by quality control tests to ensure they received the proper degree of treatment.

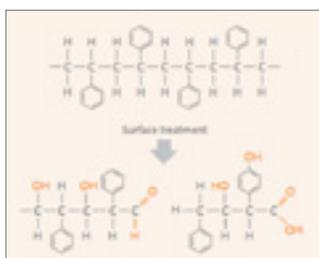


Figure 1. After treatment the hydrophobic polystyrene chain is oxidized to a more wettable, hydrophilic surface for better cell attachment.

Since these modified surfaces are not visibly different from the untreated surface, there is no easy way, short of growing cells, for customers to check the adequacy of the treatment process. As a result, many customers assume cell attachment and growth problems are caused by mistakes made during manufacturing. It is very important that the plastic culture vessel's status as the suspected cause of the problem be resolved as quickly as possible so the real cause can be identified and eliminated. Usually, the first step is to carefully read the packaging that the vessel came in to ensure that it received the standard surface treatment for cell culture. Some vessels, especially flasks, dishes, and microplates, are also available without any tissue culture treatment. These are designed for special applications where an untreated surface is required usually to prevent or reduce cell attachment. Carefully reading the packaging labels before use will prevent these mistakes.

Corning also offers several unique surface treatments to enhance cell attachment for special applications including the Corning® CellBIND® surface, Corning Primaria™, Corning PureCoat™ Amine, and PureCoat Carboxyl surfaces. For some cell types, such as primary cells or stem cells, an extracellular matrix coating may be required. These unique treatments also alter the surface charge of culture vessels, improving the attachment and growth of fastidious cell types, such as primary or transfected cell lines in reduced- or serum-free environments. Corning enhanced surfaces are suitable for research, drug discovery, and high throughput screening applications when the standard surface treatments are not enough for good cell attachment.

Next, compare the performance of the suspected vessel to the same product from a different production lot, or to similar products from another manufacturer. If a difference is found or the results are not clear, then it is time to contact the product manufacturer for assistance. Once the vessel is eliminated as the problem's cause, the customer can then continue to look elsewhere for a solution. The following examples will help identify some common problems and causes often incorrectly associated with poor surface treatment of plastic vessels and, where possible, will offer some solutions that can be used to eliminate them.

For additional information on surfaces for better cell attachment, refer to the Corning Cell Culture Surfaces Selection Guide (Corning Lit. Code CLS-C-DL-AC-006).

Problems Related to Technique

The greatest opportunity for cell culture problems occurs during the day-to-day activities of feeding and maintaining the cultures. Culture loss due to contamination is one of the most frequent and more serious of these problems. However, many other problems, while less serious and not as noticeable as contamination, still adversely affect the cultures and experiments. Often, the first sign that something is wrong occurs when the cells are microscopically examined, and uneven or unusual patterns of cell attachment or growth are observed. Identifying the problem's cause is the first and usually most difficult step, followed by finding the right solution. Sometimes transient growth problems occur and then disappear without ever identifying a cause. **NOTE:** Many of these growth problems are not readily observed during routine microscopic observation of live cultures. The occurrence and extent of these problems are best observed when sample cultures are first fixed (2.5% glutaraldehyde or 70% ethanol) and stained (1% crystal violet stain) prior to observation.

Growth Problems in Flasks, Dishes, and Plates

Poor or Slow Cell Attachment

Cell attachment problems are often serious, especially when growing cells in reduced- or serum-free medium. The attachment proteins (vitronectin and fibronectin) and other proteins found in serum containing culture media are essential for many cell lines to adhere and spread on the vessel bottom. Switching to serum-free media removes these attachment proteins forcing cell lines to start manufacturing their own attachment proteins in order for them to attach well to the vessel surface. It may take cell lines several passages before they are fully adapted. Primary cultures, especially in serum-free media, will also attach and perform better by using culture vessels pre-coated with attachment proteins or extracellular matrix. For cells that have attachment problems on traditional cell culture vessel surfaces, Corning recommends trying the Corning CellBIND surface or other enhanced treatment or coated surfaces. Pre-warming the medium used for the initial cell seeding and pre-gassing larger culture vessels, so the medium will reach its correct pH sooner will help cells attach more quickly. Pre-gassing larger vessels before seeding is highly recommended and should be done with filtered medical grade 5% CO₂/95% air mixtures.

Spotting

Clear areas, often resembling single or small clusters of viral plaques, occur along the sides or in the middle of the vessels (Figure 2). This usually results when the initial cell inoculum contains foam or bubbles. Although they appear to float on top of the medium, bubbles also push below the surface, displacing the cell-containing medium and reducing the cells available for attaching. Should the bottom of the bubbles contact the bottom of the vessel, it will prevent cells from attaching in that area. Bubbles only need to stay intact for a short period to have this effect. Bubbles that occur during refeeding of cultures, but after cells have attached, may cut off the cells from the fresh medium. If the bubbles last long enough, this will result in cell necrosis under those areas. Bubbles can usually be avoided by careful attention to mixing and pipetting techniques.

For information on preventing culture loss from contamination, refer to A Guide to Understanding and Managing Cell Culture Contamination (Corning Lit. Code CLS-AN-020).



Figure 2. Examples of spotting in flasks caused by bubbles in the medium. The flask on the bottom right shows a normal bubble-free cell monolayer. These flasks were fixed and stained to show the effects of bubbles on cell attachment and subsequent growth.

Uneven Growth

Inadequate mixing of cell inoculum with medium during addition to vessels, especially dishes, can result in uneven distribution, attachment, and growth of cells (Figure 3). Experience or simple experimentation will generally result in effective ways of mixing medium thoroughly without creating bubbles or foam.

Uneven growth can also occur as a result of the shear forces generated by medium sweeping across cell monolayers during medium changes or while moving cultures between the laminar flow hood and the incubator¹⁷. This effect is often more pronounced in serum-free cultures.

Static Electricity

Static electrical charges that build up on plastic vessels can also adversely affect cell attachment. This problem occurs more frequently when the relative humidity is very low during the winter (or year-round in some laboratory locations). Wiping the outside of vessels with a clean damp towel, increasing the room humidity, or using commercially available antistatic devices may eliminate or reduce this problem. Extra care should be used to avoid rubbing the vessels against the packaging when opening them (especially roller bottles), as this can increase the static charge.

Meniscus Rings

When too small a volume of cell inoculum is used, a pattern of heavier growth will appear along the sides of flasks or as a ring or halo in dishes. (This halo effect is often more pronounced in multiwell plates.) This pattern occurs because the meniscus along the sides of the vessel is deeper and contains more cell inoculum and medium per unit surface area than does the thinner film of inoculum toward the center of the vessel. A similar effect is observed when too little medium is used in refeeding cultures. As a general rule, 0.2 to 0.3 mL of medium should be used for every cm² of growth surface. The large meniscus, with a thin layer of media in the center, can also appear if the incubator is not properly humidified. This will not show as uneven attachment, but the increase in osmolality will impact cell growth (Table 1).

Mycoplasma

Although the problem of cell culture contamination is beyond the scope of this guide, it is important to draw attention to another potential and widespread source of mysterious cell attachment and growth problems. For more detailed information on the problem of cell culture contamination, refer to References 8, 11, 13, and 15. Due to the very high densities they can achieve in cell culture (up to 10⁸/mL), mycoplasmas (unlike other contaminants such as bacteria and fungi) cause serious adverse effects on cell cultures without clouding the medium or being observed under the microscope.



Figure 3. The cell inoculum was not adequately mixed when added to this flask. Consequently, part of the flask surface was not covered by medium when it was placed in the incubator.

For additional information on detecting and preventing mycoplasma contamination, refer to the Mycoplasma Detection Using DNA Staining protocol (Corning Lit. Code CLS-AN-025).

Mycoplasmas often grow attached to the cell membrane; as a result, a single cell may have several hundred mycoplasma on its membrane which greatly affect the ability to attach and grow. An ongoing mycoplasma screening program is an essential requirement for all cell culture labs working with cell lines^{8,10,12,13}. Without such a program, mycoplasma contamination, along with the associated problems, is likely to occur at some point. Mycoplasma screening can be done in-house or sent out for testing. It is best not to use cell-based mycoplasma tests in house. If doing in-house tests, use DNA-based detection kits.

Growth Problems in Roller Bottles

The constant movement of the medium across the surface of the bottle, as slow as it appears, can make it more difficult for cells to attach and grow in roller bottles compared to stationary vessels such as flasks and dishes. The constant motion of the medium can also lead to a more stressful cell environment than is found in stationary culture systems. Consequently, any technique-related issues that reduce the attachment ability of cells is magnified and clearly stands out⁴. **NOTE:** Many of these growth problems are not readily observed during routine microscopic observation of live cultures. The occurrence and extent of these problems are best observed when sample cultures are fixed and stained prior to analysis.

Uneven Cell Attachment and Clumping

One of the most frequently encountered problems using roller bottles is the difficulty of getting the cells to attach and form an even monolayer in the bottle. Rotating bottles at inappropriate speeds is a common cause of attachment problems. If the bottles are rotated too quickly for cells to easily attach, areas of heavy cell growth often appear as circular bands towards both ends of the bottles (Figure 4). This is because the medium flow is slightly slower at the ends than in the middle of the bottles. Rotating bottles too fast may also result in large clumps of cells. This results from the tendency of cells to form clumps since they find it easier to adhere to each other than to the surface of the roller bottle. Eventually, these clumps become large enough to attach to the bottle surface. A recommended starting speed for initiating roller bottle cultures is 0.5 to 1.0 revolutions per minute (rpm) to start. However, if cells have difficulty attaching, slower speeds (0.1 to 0.4 rpm) should be used until the cells are attached.

Cell damage during subculturing, or incomplete inactivation or removal of dissociating enzymes can also make it more difficult for cells to attach and may result in banding or clumping. The protein-based cell receptors used to initiate cell attachment become damaged by the dissociating procedures and must be replaced before the cells can reattach. Serum is commonly used to deactivate dissociating enzymes. If using serum-free media, other methods will be required. Poorly regulated incubator temperatures (temperatures that are too high or too low) will also make it more difficult for cells to evenly attach to roller bottles.

If the bottles are initially rotated too slowly, or if they slip or stop turning even for a short time during the initial cell attachment period, uneven longitudinal bands of cell growth may appear. Cleaning the rollers on the roller apparatus should alleviate slipping bottles. If necessary, rubber bands can be placed around the ends of the bottles to improve traction.

Table 1. Recommended Medium Volumes for Corning® Dishes and Flasks (mL)

Corning Plastic Culture Flasks Surface Area (cm ²)	Recommended Medium Volumes (mL)
25	5 to 7.5
75	15 to 22.5
150	30 to 45
175	35 to 52.5
225	45 to 67.5

Corning Plastic Culture Dishes Diameter (mm)	Recommended Medium Volumes (mL)
35	1.6 to 2.4
60	4.2 to 6.3
100	11 to 16.5
150	30 to 45
245 (square)	100 to 150


*Based on using 0.2 to 0.3 mL medium per cm² of growth area.



Figure 4. The roller bottle on the left was rotated too fast resulting in uneven distribution of cells to the ends of the bottle. The bottle on the right was rotated at the correct speed.



Figure 5. As a result of the roller apparatus not being level, most of the cells have attached to the bottom end of this bottle. This also can occur when bottles are stood on end for too long immediately after seeding it with cells.

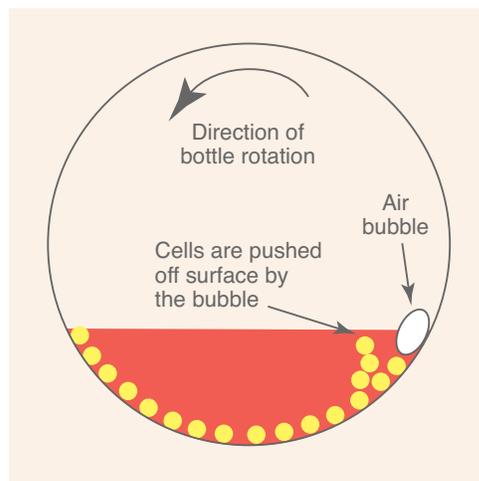

Bands of heavy growth at just one end of the bottle are often the result of the roller apparatus not being level, causing an increased amount of medium and cells at the end of the bottle that is lower (Figure 5). Furthermore, the longer it takes the cells to attach, the more time there is for them to gradually roll down the side of the bottle to the lower end before attaching. Standing a bottle on end for too long after initially seeding it with cells can have a similar effect.



Figure 6. These clear bands were caused by debris in the medium scraping away cells as the bottle was rotated.

Clear Bands

Occasionally, clear circular bands will occur on roller bottles where the cells appear to have been swept away (Figure 6). While small pieces of rolling debris or large cell clumps may cause this to occur, one of the most common causes is the short-term presence of bubbles in the initial cell inoculum. These bubbles, when in contact with the sides of the slowly rotating bottle, can act as miniature plows, scraping off the cells as they begin to attach. Avoid bubble formation by carefully pouring medium down the sides of the bottles, or by pipetting it directly into the bottom of the bottles. Cell suspensions used for inoculating roller bottles should be carefully prepared to ensure they are bubble-free.

Streaking

Condensation (essentially pure water) falling onto exposed cells can cause some unusual patterns and events. This problem usually occurs in roller bottles that have been removed from an incubator and are standing upright at cooler room temperatures while awaiting processing. Due to temperature differences, water vapor will condense on the inside of the cap. The resulting droplets may then coalesce and run down the sides of the bottle across the cells that are now only covered by a very thin film of medium. These cells will then undergo a strong osmotic shock. If they have formed a confluent monolayer, they may tear or pull apart from each other along the path the water takes, creating a visible dagger-like streak (Figure 7). Cells that have not reached confluency may round up and float off into the medium, leaving behind a long clear streak devoid of cells.



Figure 7. A clear, streak-free of cells resulting from cap condensation running down the cell sheet while the bottle was temporarily stored upright awaiting processing.

Peeling

Heavily confluent cell monolayers (especially fibroblasts) will occasionally start to peel away from the surface of the roller bottle. This also occurs in flasks, dishes, and microplates. This results from the formation of a flexible sheet of tightly interconnected cells and cell-manufactured extracellular matrix, not from surface treatment failure. Over time, mechanical stresses can develop in the cell sheet from cellular movements and contractions that may then cause the cell sheet to tear or pull away from the roller bottle (Figure 8). Physical damage from pipetting directly onto the cell sheet, tearing it with the end of the pipet, or other manipulations to the cell sheet may also initiate cell sheet peeling.

Problems Related to Incubators

Cells spend nearly all of their existence in incubators, yet these units may not always provide the stable, consistent environment cells require. Besides the obvious function of maintaining temperature, incubators used in open culture systems also control humidity, the gaseous environment around the cells, and indirectly, the pH of the culture medium. Ideally these parameters should be constant and not a source of experimental variation. Unfortunately, variation does occur and can be a major problem if not recognized and eliminated^{4,5}.

Temperature

Temperature differences within the incubator, even though small, can create problems when the differences are only a few tenths of a degree. Constant opening and closing of poorly insulated incubator doors can result in significant temperature reductions, usually localized toward the front of the incubator. Often this effect is first noticed when heavy condensation forms on vessels located near the front of the incubator. As a result of slightly cooler temperatures, these vessels may have considerably slower growth rates than their neighbors to the rear, as well as being more prone to fungal contaminants from the condensate. One solution is to set aside a separate incubator, with reduced traffic in and out, for all critical experiments, thus minimizing temperature fluctuations. Where this is not practical, those areas within the incubator that have the least temperature fluctuation (usually towards the rear) should be utilized for critical work.



Figure 8. A cell monolayer damaged by scraping with a pipet (left). Cells peeling away from the surface of the bottle (right).

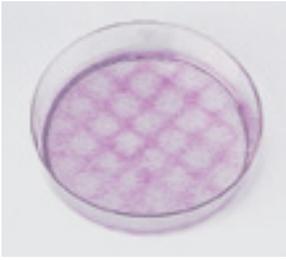


Figure 9. Growth pattern showing differential attachment and growth of WI-38 human fibroblasts. This pattern matches the pattern of holes on the shelf on which the dish was incubated.



Figure 10. Growth pattern showing differential attachment and growth of cells in a flask. Here the growth is heavier over the holes. Note also the clear area caused by bubbles.

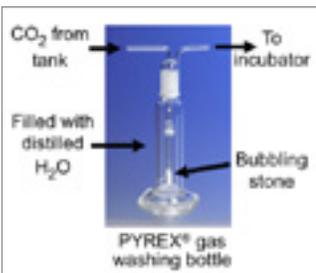


Figure 11. Humidifying incoming gases will reduce evaporation problems in incubators.

Usually temperature effects are difficult to visualize and are less likely to be recognized. Figure 9 shows a distinctive growth pattern for WI-38 cells that was caused by minor temperature fluctuations in a dish. The dish was lightly inoculated with cells, placed in an incubator on a perforated metal shelf for 48 hours, and then fixed and stained. The relatively clear areas on the plate perfectly match the perforations on the metal shelf showing the cells preference for the warmer areas directly over the metal portion of the shelf. These conditions usually occur when incubators are frequently opened, especially during the first few hours after freshly inoculated cultures are placed inside.

A similar pattern in a flask is shown in Figure 10. Here the cell growth and attachment is better over the holes in the shelving. This indicates the temperature in the incubator is a bit too high for the cells. These patterns indicate that cells are sensitive to relatively small temperature changes (less than half a degree).

Stacking vessels together can also result in vessel-to-vessel differences in temperature and growth rate. The vessel on the bottom of the stack, which is in contact with the metal shelf, warms up fastest when initially placed in the incubator. The vessel on the top is likely to cool faster, while a vessel in the middle is more insulated from any temperature fluctuations. It is very important to consider these positional effects when designing experiments where growth rates will be evaluated. To help avoid these problems, use spacers or empty “dummy” vessels to avoid direct contact with perforated shelves and if possible, do not stack the vessels.

Evaporation

Evaporation-induced changes in the osmolality of the culture medium can affect both the cell growth rate and, occasionally, the patterns of growth. Evaporation effects are easily observed in 96-well microplates where the outer peripheral wells (especially the four corner wells) often show a marked reduction in media volume over time. While tighter fitting lids can reduce this effect, it cannot be eliminated. Visible evaporation effects in other types of culture vessels, although present, are much harder to detect.

Evaporation losses depend on the type of culture vessel used (flasks, dishes, multiwell plates, etc.), their location in the incubator, and frequency of entry into the incubator. The humidification system, positioning of circulating fans, amount of insulation, and general airflow patterns will all help determine local evaporation levels. Evaporation can be minimized by keeping water reservoirs full and humidifying incoming gases (usually carbon dioxide) using a gas washing cylinder (Figure 11, Corning 31770-500EC, for example) filled with purified water in-line with any gases being fed into the incubator.

The following method can be used to create a ‘map’ for an incubator showing both the expected evaporation levels and those areas to be avoided for experiments requiring long-term incubation.

Procedure for determining evaporation levels in different sections of an incubator

1. Determine the total number of dishes that will be needed and how many shelves and dishes per shelf will be tested. At least 3 shelves (top, middle, and bottom) with 9 to 16 dishes per shelf (in a 3 × 3 or 4 × 4 matrix) are recommended for each test.
2. Consecutively number the bottoms of all dishes to be used. Then, accurately pipette the appropriate amount of medium (4 to 5 mL for a 60 mm dish) into each dish, and record the weight (without the cover in place). Water can be substituted for medium without affecting accuracy, thus eliminating some expense and the problem of contamination. Cover the dish and place in the appropriate position in the incubator. Make sure the position of each dish is recorded so the map can later be accurately constructed.
3. Incubate for the desired time period. This is usually the length of your average experiment. It is preferable during this period to continue normal use of the incubator to better reflect actual expected evaporation levels. If not, normal use should be simulated by periodically opening the incubator.



Figure 12. Vibration-induced concentric ring pattern in a 100 mm dish containing chick embryo fibroblasts that was stained 1 hour after plating.

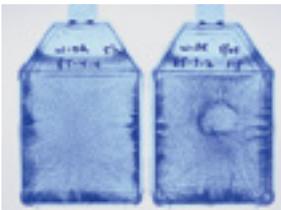


Figure 13. Flasks containing WiDr (human tumor) cells from the same incubator show varied vibration effects depending on their location in the incubator and the amount of damage the cells received during the subculturing process.



Figure 14. Wells from two 6-well plates. The well on the top shows an evenly stained cell monolayer from a plate on a level shelf. The well on the bottom shows the effect of a tilted shelf on cell growth.

- Remove the dishes in small groups from the incubator, and quickly weigh each without the covers in place. Any cover condensation represents water lost from the medium and should never be counted in the weighing process. The difference between the initial dish plus medium or water weight (in grams) and final weight, after evaporation losses, for each dish divided by the volume of water in mL initially added ($\times 100$) will very closely approximate the total evaporation loss for the dish expressed as a percent.

$$\frac{\text{Initial Weight (in grams)} - \text{Final Weight}}{\text{Volume of Water Added (in mL)}} \times 100 = \% \text{Evaporation Loss}$$

- Construct the map showing evaporation levels on the different incubator shelves. Any “hot spots” will clearly stand out; their cause can then be determined and corrected, or the area can be avoided in the future. Evaporation levels higher than 10% to 15% may have adverse effects on cultures. Wide differences in evaporation levels in different locations within the incubator will cause considerable variation in experimental results as well.

Vibration

Incubator vibration is responsible for some of the most bizarre growth patterns that occur in culture vessels. Its effect primarily occurs on cells when they are trying to initially attach to the surface following inoculation of the vessel. In dishes, vibration will push cells to the edges or the middle of the vessel, or will sometimes form concentric rings of cells. Figure 12 shows this type of pattern in a chick embryo fibroblast culture. The entire batch of 40 dishes placed in the incubator showed similar patterns. Although the vibration that caused this problem was always present in the laboratory, this pattern never occurred until the CO₂ tank supplying the incubator emptied at the same time the cultures were incubated. This loss of CO₂ raised the pH of the medium and made it more difficult for the cells to attach, allowing time for the vibration to have its effect.

Figure 13 shows typical patterns that occur in flasks because of vibration. Patterns in flasks tend to be more varied due to the irregular shape of the vessel. Changes in the frequency and amplitude of the vibration or vessel location on the incubator shelf will also affect the cell patterns. Sharpness or intensity of the resulting patterns weaken over time, especially with more motile cell lines, and are most noticeable when vessels are stained after 24 to 48 hours of incubation.

Finding the source of the vibration is difficult. It may be as simple as a loose fan motor within the incubator, but is far more likely to be caused by a more remote source. Heavy foot traffic, air handling units, compressors, laminar flow hoods, centrifuges, refrigerators, and other motorized appliances should be kept as far from incubators as possible.

Incubators should be placed on heavy, sturdy tables or benches that are not shared with any vibrating equipment. To reduce the effects of natural building vibration, position them on a floor directly over major structural supports. Spaces along well-supported outside walls, if located away from heavy traffic areas both inside and out, will often have less vibration than the central spaces. If the incubators are not fully supported and carefully leveled, then vibration effects can be enhanced. Setting up experiments at the end of the day may help eliminate some of the vibration caused by people-associated activities.

Levelness

Unusual patterns also occur when the vessels in the incubator are not level. Improper stacking of vessels or using shelves that have not been leveled prior to use often causes this. Figure 14 shows stained wells from two 6-well plates, one of which was not on a level incubator shelf. The ‘crescent moon’ effect from the lack of medium caused by the tilted shelf is clear. Shelves should be checked with a spirit or bubble level and adjustments made following the incubator manufacturer’s recommendations. It is recommended that shelves be checked periodically to make sure they are level, especially after their removal for cleaning.

Sometimes an unusual growth pattern indicates that more than a single problem is causing the pattern. Figure 15 shows a stained flask that has parallel rows of cells caused by an incubator vibration problem. However, there is also heavier cell growth on the left side of the flask indicating either that the incubator shelf was not level or the flask was improperly stacked on other flasks.



Figure 15. Vibration has caused the cells in this flask to attach in parallel rows. Moreover, the heavier cell density on the left side of this flask indicates the incubator shelf was not level.

Gases

The incubator, often considered a major source of biological contamination, can also be a source of chemical contamination. The gas mixtures (usually containing carbon dioxide to help regulate media pH) perfused through some incubators may contain toxic impurities, especially oils or other gases such as carbon monoxide, that may have been previously used in the same storage cylinder or tank. This problem is very rare in medical grade gases, but more common in the less expensive industrial grade gas mixtures. To make sure the correct gas cylinder is used, care must also be taken when installing new cylinders. Other potential chemical contaminants are the toxic, volatile residues left behind after cleaning and disinfecting incubators. Disinfectant odors should not be detectable in a freshly cleaned incubator when it is placed back into use. Most incubators have a CO₂ readout. If you are having cell growth issues, you should check the CO₂ level with a Fyrite® or similar device. For additional information on useful disinfectants see Reference 5.

Problems Related to Culture Media

As mentioned earlier, both culture media and culture vessels are prime suspects whenever mysterious cell growth or attachment problems occur without obvious causes. Unless heavily contaminated, good culture medium is not visibly different in appearance from defective culture medium. The only good way to determine medium quality is to attempt to grow cells with it; this is the basic quality control procedure used by Corning and other media manufacturers and the only good method for homemade media as well.

Depending upon both the nature and the degree of the problem, cell cultures respond to deficient or toxic media in different ways. These responses can range from minor changes in growth rate or cell attachment to the total destruction of the culture. Determining if the medium is responsible for a problem is relatively easy; simply test the suspected batch against a sample proven effective. Determining why the medium is defective is extremely difficult due to the numerous reagents and complex steps involved. Therefore, time and energy are much better spent preventing media problems than trying to find and fix them later; management by prevention is the key to successful media production. The following sections will discuss some of the common and uncommon problem areas for making and using culture media⁷.

Formulation Errors

The first step is deciding which medium is best for your applications. Good starting points are usually from scientific literature, colleagues, or from the vendor from whom the cells were sourced. Unfortunately, there is much confusion in this area from the many formulations using the same or similar acronyms. Corning offers many different formulations of 1X liquid and powdered media. Often these formulations differ from one another by omitting a single, but critical, reagent such as glutamine. It should be no surprise that confusion and mistakes are frequently made when selecting or ordering media. Read the labels on new bottles of medium very carefully, and make sure the bottle's contents are identical to previous medium and contain all the necessary components.

Formulations printed in scientific journals also contain occasional errors or intentional changes that add to the confusion³. Special care and effort must be taken to ensure the medium you make or buy is what was recommended and you know its contents.

When making media from scratch, the formulations used should be carefully checked with at least two reputable sources. The most obvious media problems arise from mistakes made during media preparation. A master formulation sheet and preparation protocol, including any special instructions or precautions, should be prepared for all media and solutions, and then used by everyone in the lab. Preparation log sheets listing all reagents used, their lot numbers, weight or volumes (both desired and delivered), pH, date, preparers, and storage conditions, are essential to insure consistency of product, reduce the potential for errors, and track down mistakes if they occur.

Reagent Quality

The next major source of problems is the medium ingredients including the water, reagents, and any special supplements such as sera and antibiotics. The highest quality water available should always be used. Special care must be taken to remove all trace metals, dissolved organics, and endotoxins. Endotoxins can have a variety of effects on cultured cells and are often found in water purified by systems that use ion-exchange resins. Purified water should not be stored for long periods before use. Water quality should be periodically checked using sensitive growth assays at clonal densities against known standards⁴.

For information on the effect of endotoxins on cell culture, refer to the Endotoxins and Cell Culture application note (Corning Lit. Code CLS-TC-305).

It is wise to purchase chemicals of the highest purity available to avoid problems associated with trace contaminants. This is especially important for serum-free media where the effects of trace contaminants are not masked by sera. To prevent any breakdown in quality, the optimum storage conditions for the chemicals must be carefully maintained once purchased. The same lot of chemicals should be used each time medium is made; when a lot is replaced, the new medium should be immediately compared against a previous batch to insure the replacement medium is satisfactory.

Buffers

After mixing, the medium is then buffered by the addition of sodium bicarbonate and adjusted to the proper pH. It is also helpful to check the osmolality at this point; mistakes in adding reagents can be uncovered as a result of finding unexpected deviations in osmolality measurements.

Many growth problems result when customers do not supply the CO₂ levels required by the bicarbonate-based buffering system of the medium they are using. This results in poor pH control and will adversely affect cell attachment and growth. For optimum buffering capacity, usually the higher the level of sodium bicarbonate, the higher the level of CO₂ is required (Table 2).

The most common system is a medium buffered with Earle's Balanced Salts containing 2.2 g/L of sodium bicarbonate. This system is designed for use in open culture vessels (dishes, microplates, or flasks with loose or vented caps) that allow gas exchange with a humidified and enriched CO₂ environment (usually 5%). The second system is a medium buffered with Hank's Balanced Salts containing only 0.35 g/L of sodium bicarbonate. This buffering system is designed for use in a sealed or gas-tight system and is not suitable for dishes and microplates.

Often, the above bicarbonate-based buffer systems are supplemented with the addition of HEPES, a widely used organic buffer. Upon exposure of the medium to fluorescent light, the use of this organic buffer can lead to additional problems. (For more information, see the section on Fluorescent Light-induced Toxicity).

Table 2. Commonly used cell culture media with the amounts of sodium bicarbonate used for buffering. Higher levels of sodium bicarbonate usually require higher levels of CO₂ added to the incubator.

Cell Culture Media	Sodium Bicarbonate Levels (g/L)	Extra CO ₂ Needed
Leibovitz's L-15 Medium	None	No
Eagle's MEM (Eagle's Minimal Essential Medium) with Hank's Balanced Salts	0.35	No
Medium 199 with Hank's Balanced Salts	0.35	No
Ham's F12	1.176	Yes
DMEM/F12	1.2 to 2.438	Yes
RPMI 1640	2.0	Yes
MEM (Eagle's Minimal Essential Medium) with Earle's Balanced Salts	2.2	Yes
McCoy's 5A	2.2	Yes
Medium 199 with Earle's Balanced Salts	2.2	Yes
MEM Medium with Earle's Balanced Salts	2.2	Yes
CMRL 1066 Medium with Earle's Balanced Salts	2.2	Yes
Iscove's Modified Dulbecco's Medium	3.024	Yes
DMEM (Dulbecco's Modified Eagle's Medium)	3.7	Yes

Filtration

Sterile filtration is usually the last step prior to the addition of any sterile supplements. There are two potential problems that may occur: first, interfering substances may be washed into the medium from the membrane; or second, valuable medium components may be lost by binding to the filter membrane. While most membranes contain usually harmless trace levels of leachable (extractable) substances, some membranes, especially some cellulose-based membranes, contain wetting agents that, at higher concentrations, may affect cell growth. These agents can be easily and safely removed from filters prior to use by running a small amount of warm high purity water through the filter unit. Some membrane materials, such as cellulose nitrate or nylon, may bind specific medium components, especially peptides and proteins used as growth or attachment factors². Testing should be performed to ensure these important supplements or factors are not being lost due to filtration.

Sera

The sera used as media additives have long been a source of problems in cell culture. These problems include the high cost of serum, its variable quality and performance, and its potential as a source of mycoplasmas, endotoxins and other contaminants. While many advances have been made in the use of serum-free and reduced-serum media, many cell culturists still use sera in their media. When practical, it is still a good idea to pre-test several lots of serum before choosing one for purchase. The most sensitive assay possible that reflects the expected use of the serum should be used.

For information on selecting the best filters for cell culture, refer to the Corning Filtration Guide (Corning Lit. Code CLS-FIL-004).

Fluorescent Light-induced Toxicity

The deleterious effect of fluorescent light on culture media may be the single most overlooked source of chemically induced cytotoxicity. It is very important to store media and cells growing in culture vessels in the dark away from sources of fluorescent light that will interact with light sensitive media components (riboflavin, tryptophan, and HEPES). These interactions result in the production of hydrogen peroxide and free radicals that are directly toxic to cells. This well-documented problem is often ignored when there are cell growth issues^{18,19}. Since the toxic effects of improperly stored media slowly increase with time, this problem is particularly difficult to identify. Besides direct cytotoxicity, other light-induced damaging effects include genetic damage (increase in mutation rates and chromosomal aberrations).

Troubleshooting Suggestions

While it is common to experience occasional minor problems when growing cells, more frequent or reoccurring cell culture problems may seriously interfere with a research laboratory and are very costly. Because of the complex nature of cell culture, identifying the causes of these cell culture problems is often a difficult and time-consuming task. The following information will provide a framework for identifying common cell culture problems and finding solutions to them⁵.

1. The first step is to begin the process of identifying and defining the exact nature and extent of the cell culture problem(s). It is usually advisable to enlist the assistance of everyone in the laboratory or research group whose knowledge or experience might contribute any information related to this problem or who might be affected by the problem. A problem well-defined is a problem half solved.
 - a. Try to place the problem into a single category, such as poor cell growth or attachment, contamination, or one of the other areas that are detailed below in this brochure.
 - b. Determine the extent of the problem. Does the problem only affect a single cell line or is it occurring with other cell lines as well? Does the problem only affect one worker's cultures or is it affecting the cultures of others in the laboratory?
 - c. It often helps to break up complex problems into smaller pieces that can be handled and understood more easily.
2. Collect, list, and organize all the known facts surrounding the problem. Be specific, and try to find possible cause and effect relationships. Do not overlook the obvious. Do not try to solve the problem at this point in the information gathering process.
 - a. What is the temporal nature of the problem? What has changed or is new in the laboratory? Is the problem on-going or is it sporadic? Are its occurrences related to a time of day, week, or month?
 - b. Is the problem associated with the use of a piece of equipment, such as an incubator, laminar flow hood, centrifuge, or water bath? Organize all known facts surrounding the problem in the order of their probability. Be specific, look for cause and effect relationships, and then discard all facts that clearly do not apply and work with the rest. Don't overlook the obvious.
3. Once there is a clear understanding of the nature and extent of the problem, begin looking for causes. Resist the urge to fix the problem by changing things before the source of the problem is clear. This often masks the original problem and makes the problem more difficult to solve.
 - a. Identify and list any changes in the laboratory that have occurred either during or immediately preceding the start of this problem. Especially look for changes in the medium (and its serum and additives) and dissociating solutions used to grow and subculture the cells, as well as their incubator. But do not overlook culture vessels, other supplies, equipment, personnel, and procedures that can be directly linked with the cultures that are experiencing the problem(s). Good record keeping is very helpful for this step.

- b. Don't forget to consider that culture problems may be related to the presence of cryptic microbial contaminants such as mycoplasma, or chemical contaminants that are difficult to detect except through the adverse effects they cause. Mycoplasma contaminates at least 15% to 25% of all the cultures currently in use in the United States.
 - c. For additional ideas, review the examples of common cell culture problems that are discussed in the technical information sections following these general problem-solving suggestions.
4. Brainstorm and make a list of all the possible causes that can be identified. Then from this list select the best possibilities and begin evaluating them, ideally by running controlled experiments (whenever possible), until the actual cause is uncovered. Be creative. This may require some testing and experimentation.
 5. Eliminate possible causes by running a controlled experiment. For example, if it could be the incubator, place cultures in separate incubators. If it could be cells, thaw a vial of fresh cells. The same approach can be used with media, water baths, etc.
 6. Fix the problem and make the necessary changes so that the chance of a future re-occurrence is eliminated or greatly reduced.

For additional assistance in this area, visit www.corning.com/lifesciences, or contact Corning Scientific Support at 1.800.492.1110; outside the United States, call +1.978.442.2200; or email ScientificSupport@corning.com.

References

1. Amstein CF, Hartman PA. Adaptation of Plastic Surfaces for Tissue Culture by Glow Discharge. *J. Clin. Microbiol.* 2:46-54, 1975.
2. Brock TD. *Membrane Filtration: A User's Guide and Reference Manual*. Madison, Wisconsin: Science Tech, 1983.
3. Burke CN, Croxall G. Variation in Composition of Media and Reagents Used in the Preparation of Cell Cultures from Human and Other Animal Tissues: Dulbecco's, Earle's, and Hank's Balanced Salt Solutions. *In Vitro* 19: 693-698, 1983.
4. Freshney RI. *Culture of Animal Cells: A Manual of Basic Technique*. New York: Allan R. Liss, Inc. Third Edition, 1994.
5. Geraghty RJ, Capes-Davis A, Davis JM, Downward J, Freshney R, Knezevic I, Lovell-Badge R, Masters JRW, Meredith J, Stacey GN, Thraves P, and Vias M. Guidelines for the Use of Cell Lines in Biomedical Research. *British Journal of Cancer* (2014) 111, 1021-1046. (Available online).
6. Hudis M. Plasma Treatment of Solid Materials. In Hollahan JR, Bell AT, Ed. *Techniques and Applications of Plasma Chemistry*. New York: John Wiley and Sons, 113-147, 1974.
7. Jakoby WB and Pastan JH, Ed. *Methods in Enzymology: Cell Culture*. (Especially Chapters 1, 2, 5, 7, and 10) Vol. 58, New York: Academic Press, 1979.
8. Lincoln CK and Gabridge MG. Cell Culture Contamination: Sources, Consequences, Prevention and Elimination. In *Animal Cell Culture Methods*, edited by JP Mather and D Barnes, *Methods in Cell Biology*, Volume 57, Chapter 4, Academic Press, San Diego, 49-65, 1998.
9. MacMichael GJ. The Adverse Effects of UV and Short Wavelength Visible Radiation on Tissue Culture. *Am. Biotech. Lab*, July/August, 1986.
10. Masover GK and Buck DW. Detecting Mycoplasmas in Cell Cultures. *MD & DI*: 43-98, 1983.
11. McGarrity GJ. Spread and Control of Mycoplasma Infection of Cell Cultures. *In Vitro* 12:643-647, 1976.
12. McGarrity GJ, Sarama, J; and Vanaman, V. Cell Culture Techniques. (Excellent overview of mycoplasma contamination) *ASM News* 51:170-183, 1985.
13. McGarrity GJ. Detection of Mycoplasma Infection of Cell Cultures. In K. Maramorosch; Ed. *Advances in Cell Culture*. Vol. 2, New York, Academic Press, 99-132, 1982.
14. Ramsey WS, Herti W, Nowlan ED, and Binkowski NJ. Surface Treatments and Cell Attachment. *In Vitro* 20: 802-808, 1984.
15. Rottem S and Barile MF. Beware of Mycoplasmas. *Trends Biotechnol* 11:143-150, 1993.
16. A Guide to Understanding and Managing Cell Culture Contamination (Corning Lit. Code CLS-AN-020) available at www.corning.com/lifesciences.
17. Tchao R. Fluid Shear Force and Turbulence-Induced Cell Death in Plastic Tissue Culture Flasks. *In Vitro Toxicol* 9:93-100, 1996.
18. Wang RJ. Effect of Room Fluorescent Light on the Deterioration of Tissue Culture Medium. *In Vitro* 12: 19-22, 1976.
19. Wang RJ and Nixon BT. Identification of Hydrogen Peroxide as a Photoproduct Toxic to Human Cells in Tissue Culture Medium Irradiated with "Daylight" Fluorescent Light. *In Vitro* 14:715-722, 1978.
20. Waymouth C. Osmolality of Mammalian Blood and of Media for Culture of Mammalian Cells. *In Vitro* 6: 109-127, 1970.

Ordering Information

Corning® Cell Culture Flasks

Cat. No.	Surface Area (cm ²)	Flask Style	Neck Style	Cap Style	Qty/Pk	Qty/Cs
430639	25	Rectangular	Canted	Vent	20	200
430641U	75	U-shaped	Canted	Vent	5	100
430825	150	U-shaped	Canted	Vent	5	50
431080	175	U-shaped	Canted	Vent	5	50
431082	225	Rectangular	Canted	Vent	5	25
10020	1,720	Corning HYPERFlask®	-	-	4	4
10030	1,720	Corning HYPERFlask®	-	-	1	4

Costar® Multiple Well Plates, Tissue Culture-treated

Cat. No.	Description	Plate Type	Qty/Pk	Qty/Cs
3516	6-well	Clear	1	50
3513	12-well	Clear	1	50
3526	24-well	Clear	1	50
3548	48-well	Clear	1	100
3596	96-well	Clear	1	50
3595	96-well	Clear*	1	50
3599	96-well	Clear	1	100
3916	96-well	Black	20	100
3917	96-well	White	20	100
3603	96-well	Black/Clear-bottom	1	48
3610	96-well	White/Clear-bottom	1	48

*Low evaporation lid

Corning Dishes, Tissue Culture-treated

Cat. No.	Dish Style* (mm)	Approx. Height (mm)	Growth Area (cm ²)	Qty/Pk	Qty/Cs
430165	35	10	8	20	500
430166	60	15	21	20	500
430167	100	20	55	20	500
430599	150	25	148	5	60
431110	245	25	500	4	16

*Dish Style (mm) 35 mm = 34.4 mm; 60 mm = 52.1 mm; 100 mm = 83.9 mm; 150 mm = 139.1 mm. Square dishes have interior bottom dimensions 224 x 224 mm.

Corning Cell Scrapers

Cat. No.	Description	Blade Length (cm)	Handle Length (cm)	Qty/Pk	Qty/Cs
3008	Cell lifter	1.9	18	1	100
3010	Scraper, small	1.8	25	1	100
3011	Scraper, large	3	39	1	100

Corning Cell Strainers

Cat. No.	Description	Qty/Pk	Qty/Cs
431750	40 µm, blue	1	50
431751	70 µm, white	1	50
431752	100 µm, yellow	1	50

Corning Cryogenic Vials

Cat. No.	Capacity (mL)	Style	Self-standing	Qty/Pk	Qty/Cs
430659	2	Round bottom, internal thread	Yes	50	500
430488	2	Round bottom, external thread	Yes	50	500

Falcon® Cell Culture Flasks, Tissue Culture-treated

Cat. No.	Surface Area (cm ²)	Flask Style	Neck Style	Cap Style	Qty/Pk	Qty/Cs
353018	25	Rectangular	Canted	Vent	20	100
353136	75	Rectangular	Canted	Vent	5	60
355001	150	Rectangular	Canted	Vent	5	40
353112	175	Rectangular	Straight	Vent	5	40
353138	225	Rectangular	Canted	Vent	5	30
353143	525	3-Layer	-	Vent	2	12
353144	875	5-Layer	-	Vent	1	8

Falcon Cell Culture Plates, Tissue Culture-treated

Cat. No.	Description	Plate Type	Qty/Pk	Qty/Cs
353046	6-well	Clear	1	50
353043	12-well	Clear	1	50
353047	24-well	Clear	1	50
353078	48-well	Clear	1	50
353072	96-well	Clear	1	50

Falcon Cell Culture Dishes

Cat. No.	Dish Style (mm)	Approx. Height (mm)	Growth Area (cm ²)	Qty/Pk	Qty/Cs
353001	35	10	11.78	20	500
353002	60	15	21.29	20	500
353003	100	20	58.95	20	200
353025	150	25	156.36	10	100

Falcon Cell Strainers

Cat. No.	Description	Qty/Pk	Qty/Cs
352340	40 µm, blue	1	50
352350	70 µm, white	1	50
352360	100 µm, yellow	1	50

Falcon Cell Scrapers

Cat. No.	Description	Qty/Pk	Qty/Cs
353085	1.8 cm TPE blade with 18 cm polystyrene handle	1	100
353086	1.8 cm TPE blade with 25 cm polystyrene handle	1	100
353089	3.0 cm TPE blade with 25 cm polystyrene handle	1	100
353087	3.0 cm TPE blade with 40 cm polystyrene handle	1	100

Corning® CoolCell® Containers

Cat. No.	Description	Capacity (Vials)	Exposed Vial Tops	Qty/Pk	Qty/Cs
432000	CoolCell, purple	12	No	1	1
432001	CoolCell LX, purple	12	Yes	1	1
432002	CoolCell LX, green	12	Yes	1	1
432003	CoolCell LX, orange	12	Yes	1	1
432004	CoolCell LX, pink	12	Yes	1	1
432138	CoolCell LX, 4 colors (purple, green, orange, pink)	12	Yes	--	4
432005	CoolCell 5 mL LX, purple	12	Yes	1	1
432006	CoolCell FTS30, purple	30	Yes	1	1
432007	CoolCell FTS30, orange	30	Yes	1	1
432008	CoolCell FTS30, green	30	Yes	1	1
432009	CoolCell FTS30, pink	30	Yes	1	1
432010	CoolCell SV2	12	Yes	1	1
432011	CoolCell SV10	6	Yes	1	1

Corning Vacuum Filtration Systems

Cat. No.	Membrane	Funnel/Bottle Volume (mL)	Pore Size (µm)	Qty/Cs
431153	PES	150/150	0.22	12
431096	PES	250/250	0.22	12
431097	PES	500/500	0.22	12
431098	PES	1,000/1,000	0.22	12

Corning Bottle Top Vacuum Filtration Systems

Cat. No.	Membrane	Volume (mL)	Neck Size (mm)	Pore Size (µm)	Qty/Cs
431161	PES	150	45	0.22	48
431118	PES	500	45	0.22	12
431174	PES	1,000	45	0.22	12

Corning Syringe Filters

Cat. No.	Diameter (mm)	Pore Size (µm)	Housing Material	Membrane Material	Sterile	Inlet/Outlet	Packaging	Qty/Cs
431219	28	0.2	SFCA	AC	Yes	LL/LS	Indiv.	49
431222	25	0.2	RC	PP	Yes	LL/LS	Indiv.	50
431224	25	0.2	NY	PP	Yes	LL/LS	Indiv.	50
431229	20	0.2	PES	AC	Yes	LL/LS	Indiv.	50

Reagents

Cat. No.	Description	Size	Qty/Cs
25-950-CQC	Dimethyl Sulfoxide	250 mL	1
354253	Cell Recovery Solution	100 mL	1
354216	Calcein AM fluorescent dye, 500 µg (10 x 50 µg)	500 mg	1
354217	Calcein AM fluorescent dye	1 mg	1
354218	DiIC12(3) fluorescent dye	100 mg	1
25-900-CI	Trypan Blue Solution, 0.4% (w/v) in PBS, pH 7.5 ± 0.5	100 mL	1
354235	Dispase	100 mL	1

Sera

Cat. No.	Description	Size	Qty/Cs
35-010-CV	Fetal Bovine Serum, regular	500 mL	1
35-015-CV	Fetal Bovine Serum, premium	500 mL	1

Corning Media

Cat. No.	Description	Size	Qty/Cs
10-013-CV	DMEM [+], 4.5 g/L glucose, L-glutamine, sodium pyruvate	500 mL	6
10-017-CV	DMEM [+], 4.5 g/L glucose, L-glutamine [-], sodium pyruvate	500 mL	6
15-013-CV	DMEM [+], 4.5 g/L glucose, sodium pyruvate [-], L-glutamine	500 mL	6
15-090-CV	DMEM/F12 50:50 mix [-], L-glutamine	500 mL	6
10-092-CV	DMEM/Ham's F-12 50/50 mix [+], L-glutamine, 15 mM HEPES	500 mL	6
16-405-CV	DMEM/Ham's F-12 50/50 mix [+], L-glutamine [-], phenol red	500 mL	6
15-016-CV	Iscove's Modification of DMEM [+], 25 mM HEPES, [-], β-thioglycerol, β-mercaptoethanol, L-glutamine	500 mL	6
10-016-CV	Iscove's Modification of DMEM [+], L-glutamine, 25 mM HEPES [-], β-thioglycerol, β-mercaptoethanol	500 mL	6
10-080-CV	Ham's F-12 Medium [+], L-glutamine	500 mL	6
10-060-CV	Medium 199 (Mod.) [+], Earle's salts, L-glutamine	500 mL	6
10-022-CV	MEM alpha medium [+], Earle's salts, ribonucleosides, deoxyribonucleosides, L-glutamine	500 mL	6
15-012-CV	MEM alpha medium [+], Earle's salts [-], ribonucleosides, deoxyribonucleosides, L-glutamine	500 mL	6
10-010-CV	MEM [+], Earle's salts, L-glutamine	500 mL	6
10-009-CV	MEM [+], 1.5 g/L sodium bicarbonate, NEAA, L-glutamine, sodium pyruvate	500 mL	6
15-010-CV	MEM [+], Earle's salts [-], L-glutamine	500 mL	6
17-305-CV	MEM [+], Earle's salts [-], L-glutamine, phenol red	500 mL	6
10-040-CV	RPMI 1640 [+], L-glutamine	500 mL	6
15-040-CV	RPMI 1640 [-], L-glutamine	500 mL	6
17-105-CV	RPMI 1640 [-], L-glutamine, phenol red	500 mL	6
10-041-CV	RPMI 1640 [+], L-glutamine, 25 mM HEPES	500 mL	6
15-041-CV	RPMI 1640 [+], 25 mM HEPES [-], L-glutamine	500 mL	6

Corning Buffers

Cat. No.	Description	Size	Qty/Cs
21-040-CV	PBS (Phosphate-Buffered Saline), 1X without calcium and magnesium	500 mL	6
21-020-CV	HBSS (Hanks' Balanced Salt Solution), 1X with calcium and magnesium	500 mL	6
21-023-CV	HBSS, 1X with calcium and magnesium without phenol red	500 mL	6
21-021-CV	HBSS, 1X without calcium and magnesium	500 mL	6
21-022-CV	HBSS, 1X without calcium, magnesium, and phenol red	500 mL	6
21-030-CV	DPBS (Dulbecco's Phosphate-Buffered Saline), 1X with calcium and magnesium	500 mL	6
21-031-CV	DPBS, 1X without calcium and magnesium	500 mL	6
25-015-CI	Corning glutagro™ supplement	100 mL	1
25-021-CI	Trace Elements A, 1,000X	100 mL	1
25-022-CI	Trace Elements B, 1,000X	100 mL	1
25-023-CI	Trace Elements C, 1,000X	100 mL	1
25-030-CI	MEM amino acids, 50X	100 mL	6
25-025-CI	MEM Nonessential amino acids, 100X	100 mL	6
25-800-CR	ITS (Insulin-Transferrin-Selenium)	10 mL	1

For more specific information on claims, visit the Certificates page at www.corning.com/lifesciences.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

CORNING

Corning Incorporated Life Sciences

836 North St.
Building 300, Suite 3401
Tewksbury, MA 01876
t 800.492.1110
t 978.442.2200
f 978.442.2476

www.corning.com/lifesciences

ASIA/PACIFIC

Australia/New Zealand
t 61 427286832

Chinese Mainland
t 86 21 3338 4338
f 86 21 3338 4300

India
t 91 124 4604000
f 91 124 4604099

Japan

t 81 3-3586 1996
f 81 3-3586 1291

Korea

t 82 2-796-9500
f 82 2-796-9300

Singapore

t 65 6572-9740
f 65 6735-2913

Taiwan

t 886 2-2716-0338
f 886 2-2516-7500

EUROPE

CSEurope@corning.com

France

t 0800 916 882
f 0800 918 636

Germany

t 0800 101 1153
f 0800 101 2427

The Netherlands

t 020 655 79 28
f 020 659 76 73

United Kingdom

t 0800 376 8660
f 0800 279 1117

All Other European Countries

t +31 (0) 206 59 60 51
f +31 (0) 206 59 76 73

LATIN AMERICA

grupoLA@corning.com

Brazil

t 55 (11) 3089-7400

Mexico

t (52-81) 8158-8400

For additional product or technical information, visit www.corning.com/lifesciences or call 800.492.1110. Outside the United States, call +1.978.442.2200 or contact your local Corning sales office.